



PROGRAMMABLE DELAY AND EXPOSURE INTERVALS

NEUTRAL DENSITY CAPABILITY

CONTINUOUS OR LOOPED RUN

SERIAL, USB AND TTL INTERFACES

FULL SMARTSHUTTER™ SUPPORT



(Shown with optional IQ35-SA and stand)

LAMBDA SC SMARTSHUTTER™ CONTROLLER

The **Lambda SC** is a full-featured microprocessor controlled shutter driver for the **SmartShutter™**. It supports the same shutter command set used for our other Lambda controllers and offers enhanced performance through additional modes of operation. The additional operating modes allow programmable control over the exposure time, and the delay between trigger and shutter opening.

In the Lambda 10 family of filter wheel controllers, shutter timing is controlled directly by the time of occurrence of commands or TTL input. In most cases, the system software used with these controllers has provisions for controlling the timing of the shutter. The **Lambda SC** shutter controller may be used in applications that would benefit from enhanced control options built into the controller. The controller can be operated remotely from a computer through the USB port, or opened and closed using a logic level input.

In the free running mode, the shutter will open and close repeatedly. The time between closing and reopening can either be directly timed, or set as an interval for the entire cycle. Delay and exposure intervals can be set by the user for up

to 5 hours with millisecond resolution. This mode can be set to continuously run or loop for a specific number of cycles. The mode is initiated from power up, or a trigger pulse, and can be interrupted by using a stop command.

Because our patented **SmartShutter** incorporates a microprocessor controlled motor/wiper design, the trajectory of each move can be controlled and the motion of the blade optimized for speed or for smoothness. When speed is the most important condition, the user can select the fast mode. If the modest amount of vibration in the fast mode is objectionable and speed is not important, the "soft" mode can be selected, moving the blade slower through the overall travel.

The neutral density mode results in a partial opening of the **SmartShutter**. Any one of the 144 steps from a fully closed to fully open can be selected. Using this variable aperture opening along with a liquid light guide acts to spatially homogenize the light to produce a uniform spatial illumination that's independent of the geometry of the input illumination.

CONTROLLER

LB-SC

Lambda SC control unit, USB and power cables and manual SMARTSHUTTER™

IQ25-SA*

25mm **SmartShutter** with stand-alone housing

IQ25-LS

25mm **SmartShutter** for mounting in Lambda LS

IQ25-DG

25mm **SmartShutter** for mounting in DG-4/DG-5

IQ35-SA

35mm **SmartShutter** with stand-alone housing

IQ50-SA

50mm **SmartShutter** with stand-alone housing

* Where vignetting may be an issue, we recommend the 35mm shutter.

Mounting adapters for Nikon, Zeiss, Leica and Olympus microscopes are available.

Patent No. 7,253,575

RoHS Compliant



SUTTER INSTRUMENT

One Digital Drive • Novato • CA 94949 • Phone 415.883.0128
Fax 415.883.0572 • Web www.sutter.com • Email info@sutter.com